



Air Master Systems Standard Paint Booth

- Standard features:
 - Vapor Proof light (bulb included)
 - Exhaust Fan
 - Filter(s) (MARV 8 rating)
 - Light and Blower switches (optional)
 - Unit is Pre-wired to junction box when equipped with switches

Product Specifications:

- The AMS Paint Booth is a light duty fume hood intended for use with experiments involving low volumes of mildly irritating vapors, light smoke or paint.
- The AMS Paint Booth is not intended to be used as a ductless fume hood. In addition, this fume hood may not be used for the following applications:
 - Explosion Proof
 - Radioisotope
 - Perchloric Acid
 - High Temperature Enclosure
 - Acid Digestion Enclosure

Operation, Care, and Maintenance:

- The blower must be on and operating whenever the hood is in use.
- The AMS Paint Booth is not to be used as a storage cabinet.
- Only materials directly involved in the experiment should be placed inside the hood chamber.
- Keep the light fixture clean.
- Maintenance procedures consist primarily of clean up and replacement of worn, damaged, or non-functioning filters, and checking the exhaust capacity periodically.
- Spills should be flushed immediately using neutralizing compounds as required.